Anuário de
Químicos

asecos®

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asecos GmbH

Your personal documentation to the asecos safety cabinet

Dear Customer,

You have made a decisive investment in safety for your company by purchasing
this asecos safety cabinet.

You now own an innovative product made of top-quality materials guaranteeing
the highest quality standards.

asecos safety cabinets have complete authorisation documents. We archive the
authorisation documents for every individual cabinet, keeping them ready for you
should you ever need them (e.g. for a works inspection or similar).
Simply request them using this form.

Tear off or copy that page and return to us by fax with your address and
serialnumber of the cabinet on it.

Yours sincerely

asecos GmbH

Firma	/ Company
Abteilung	/ Department
Name	/ Name
Straße	/ Street
PLZ	/ Postal Code
Ort	/ City
Schrank Seriennummer(n)	/ Cabinet Serial-Number(s)

OPERATING INSTRUCTIONS

Dear Customer,

Thank you for purchasing this cabinet for the storage of hazardous materials. In doing so you have made a decisive investment in safety for your company.

Using our cabinets for the storage of hazardous materials makes the storage of hazardous materials at the workplace safe and easy.

Please read through these operating instructions very carefully. They explain in detail the advantages and ease of using our cabinets for the storage of hazardous materials.

This will simplify the everyday handling of hazardous materials.

Many thanks.

Your ascos Team

Table of contents of these operating instructions

Page	Section	Content
13	1	Range of models
13	2	Data
13	3	Intended purpose
13	4	General safety information/warranty
13	5	Technical data
14	6	Transport
14-15	7	Installation, conditions of installation, functional check
15	8	Closing
15	9	Door open arrest system with thermal release system
15	10	Locking
15-16	11	Interior fittings
16	12	Earthing
16	13	Ventilation
17	14	Storage volume/combined storage/provision
17	15	Marking
17	16	Opening the cabinet after a fire
17	17	Disposal
17	18	Safety check

1. Range of models

Refer to Table in Annex 1 at the end of the Operating Instructions.

2. General Data

The cabinet data is contained in the board book in the cabinet.

3. Intended purpose

Models Fxx.xxx.xxx, VBF.xxx.xxx(+)
are intended for storing flammable liquids in working areas in accordance with applicable national regulations.

Models TRG.205.xxx, G90.xxx.xxx
are intended for the provision and emptying of pressurised gas cylinders in buildings in accordance with applicable national regulations.

4. General safety information / warranty

- Basically all type of work on the electrical installation (VBFA196.xxx) has to be made only when electric is off. Therefore refer to the accident prevention prescriptions and the regulations of your provider.
- Please observe the laws and regulations applicable to the handling of hazardous materials and the information in these operating instructions.
- Only use the cabinets and accessories if they are in a proper condition.
- Only use the cabinet after instruction; access should be denied to unauthorised persons.
- The area of radius of the doors should always be kept free.
- The doors/drawers should be kept closed.
- Please observe the maximum limits to quantities stored, loading, etc.
- The sump of safety cabinets which comply with EN 14470 part 1 must have a minimum collecting volume of 10% of all packs stored in the cabinet, or at least 110% of the volume of the largest individual pack, depending on which volume is largest.
- Packs with a volume larger than the volume of the sump should not be placed in the cabinet.
- Any leaking hazardous materials should be mopped up and removed immediately.
- Containers with aggressive chemicals (acids and alkalis) must be stored in special acid or alkali cabinets, or in acid or alkali compartments.
- Before storing check the resistance of the cabinet surface to the chemical.
- The storage of corrosive fluids or gases may have an effect on the function of the shut-off devices for the ventilation air.
- Conditions of installation may require the cabinet to be bolted to the building.
- Follow the instructions of the Technical Supervisory Service.
- Ensure that the requisite safety checks are made by authorised skilled staff only.
- Incorrect function, damage, corrosive damage arising for incorrect transport and installation due to the great weight and the bulkiness of the cabinets can be avoided by trained and authorised staff.
- The cabinet must be given a safety check before use.
- **Warranty for this product is agreed between you (the customer) and your specialised dealer (the vendor). Independent of this, the manufacturer assumes a warranty of 24 months starting from date of purchase for the products listed in the Operating Instructions. As safety equipment, all models are subject to an annual obligatory inspection by skilled personnel authorised by the manufacturer. If the inspection is not performed on an annual basis or not by skilled personnel authorised by the manufacturer, the warranty claim of the customer shall be forfeited vis-à-vis the manufacturer.**

5. Technical data

Refer to Table in Annex 1 at the end of these Operating Instructions.

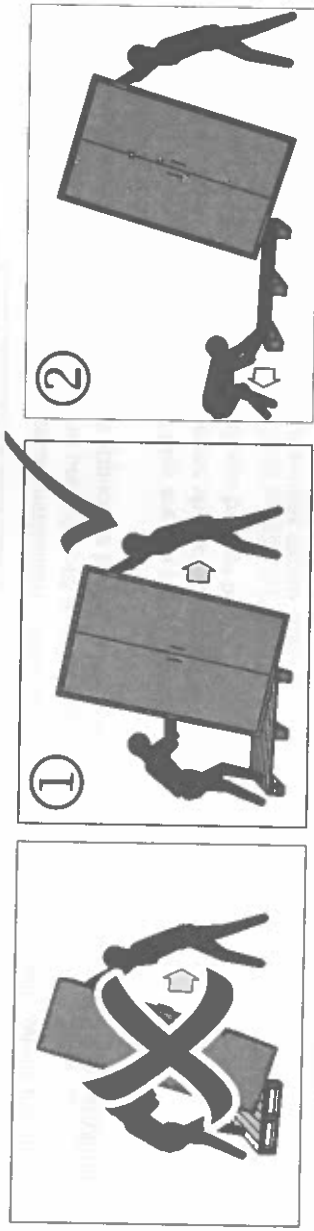
6. Transport

Take the heavy weight of the cabinets into consideration ~ refer also to the Table in Annex 1. The transport work should therefore usually be performed with a minimum of 2 people.

Transport the cabinet in an upright position using a fork lift truck; it should be secured against slipping until it reaches its final place of use.

Warning: The spacers in the door joints may only be removed at the site of installation.

The cabinet may not be laid on doors or drawers.
If the cabinet needs to be moved to its side during transport, this may only happen smoothly without jerking.


N.B.: Due to its heavy weight, the cabinet is difficult to handle. Incorrect transport may lead to concealed damage to the fire protection insulation.

We can only guarantee the requisite quality if the cabinet is transported by our specially trained staff to the point of use.

Attention: Written notification of any damage to the cabinet must be given immediately!

7. Installation/commissioning/functional check

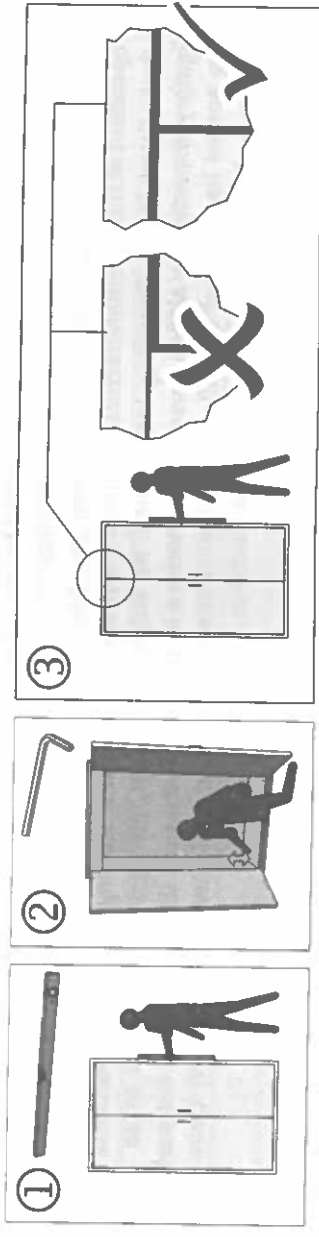The cabinet must be erected at its final site in such a way that the doors close automatically from any position and interlock correctly (with the exception of the TRG-, G90 and VBF.60 models, as well as F30.60-models with drawers). This must still be the case even if the lock has already interlocked! **When closing, the doors may not rub on the fire protection seals in the door fold!**

Metal spacers must be used to adjust the cabinet on uneven ground.

Before use, the cabinet must be given a safety inspection. Please refer to the applicable national regulations on the conditions of installation for cabinets for storage of hazardous materials.

Models F60.xxx.xxx, VBF(T/A)196.xxx(+), VBF129.xxx(+), VBF1xxx, TRG.205.xxx, G90.xxx.xxx

Cabinets of this range have 4 adjustment feet on the bottom of the cabinet for adjustment purposes (max. 10mm height adjustable). A suitable Allan key is enclosed!

Models F30.60.xxx, VBF.60.xxx

The use of a securing bracket (Art. No. EP.ML.6323) is essential to prevent the cabinet tipping over when full to maximum and when the drawers are fully extended.

Models V8F.221.xxx

The drum cabinet is supplied on wooden transport blocks. These must be removed during installation. The **spacers may not be removed**, however. First screw the enclosed brackets to the side of the cabinet. The wooden transport blocks can now be removed after lifting sideways using a fork-lift truck.

Warning: Then slowly lower the cabinet to avoid damage.

8. Closing

In the case of fire, the doors and drawers of all models (with the exception of the TRG- and G90- models) close automatically. (Cabinets of model series VBF(T) 129/196.xxx+ close automatically after 60 seconds).

If the contents of the cabinet are not being accessed, the operator/user must ensure that all doors and drawers are kept closed.

Models VBFA196.xxx

The doors of these models open and close fully automatically. The integrated sensor technology opens the door/doors automatically after personnel ID recognition. (The ID recognition is made after approximately 2 seconds). After leaving the recognition fields, the door/doors close(s) automatically after approximately 30 seconds. There is no danger of injury in the process of closure because the cabinet does not close by motor force.

Meaning of the LED displays:

ed – person is recognised, cabinet opens

Technical data:

power connection 230V, 50Hz with 5m power cable including mains plug.

9. Door open arrest system with thermal release system in the event of fire

In the case of doors with automatic closure, it is possible to fix the door open by means of a door stop system. In the case of fire, the doors close automatically due to an integrated thermal release system. (Cabinets of model series BF(T) 129/196.xxx+ close automatically after 60 seconds, door open arrest system is standard here).

0. Locking

All cabinets can be locked using a safety cylinder lock. The models of the type series VBF(T/A) 196.xxx(+), VBF129.xxx(+), BFT.xxx(+), and TRG.205.xxx and G90.xxx.xxx are equipped with a replaceable profile cylinder (possible to integrate in master key system).

B.: It should generally be remembered for all models that the cabinets do not have an emergency unlocking device, i.e. any persons locked in the cabinet cannot get out alone.

1. Interior fittings (1)

carrying force and catch volume of shelves and collecting sumps are provided in Annex 1 – Technical Data.

.1 Shelves

These shelves are already fitted to the cabinet in the factory and each secured by 4 screws. These screws must first be loosened before adjusting the height.

Warning: after adjusting the shelves, secure them again against tipping up with all 4 screws.

11. Interior fittings (2)

11.2 Bottom collecting sump/perforated metal plate insert

The bottom collecting sump, or drawers serves to collect any leaking fluids and may only be used in combination with a suitable perforated metal plate insert. In the case of a leak, the liquid in the floor collecting tray must be mopped up with suitable means, the selection of which is up to the user.

11.3 Drawers/2nd storage level

All drawers have been designed in such a way that they are permanently drawn in the cabinet or automatically drawn into the cabinet in the case of fire, making it possible to close the cabinet doors. Drawers may only be adjusted in height by authorised ascos staff as the safe drawing back of the drawers in the cabinet in case of a fire has to be guaranteed!

11.4 TRG and G90-models

- Assembly rails
The horizontal assembly rails serve to accommodate the pressurised gas fittings. The assembly rails can be height-adjusted as follows:
 - Slightly loosen screws, reposition the assembly rail and retighten screws.
- Bottle holder
is permanently attached to the rear wall of the housing and has a belt for fixing of the gas bottle(s).
- Fold away ramp
has been fitted in the factory and automatically stops in closed position. To use, simply pull the flap down and push down completely. Return to cabinet after use.
- Pipe conduit

The head part of the cabinet has fittings for a large number of pipe conduits.

Drill through the head part from the top for pipe conduits in one of the preset positions with exactly the diameter of the pipe to be fitted. The pipe can now be introduced without any further insulation measures.

- Ermeto screwed connection

The preparation for Ermeto screwed connections is made by the factory on the ceiling of the cabinet. Commercially available Ermeto screwed connections can then be fitted through the R3/8" internal thread.

12. Earthing

All models can be connected via an earthing connection in the head part or on the rear wall to an equipotential bonding.

Earthing during the decanting

Cabinets in which refilling work can be performed should additionally be equipped with an earthing terminal (upon request and with extra charge) which must be connected with the refilling containers.

Attention! In this case it is imperative to connect the external earthing connection at the safety storage cabinet to an equipotential bonding.

13. Ventilation

All models can be connected to a technical exhaust air system via the extraction air outlet (DN75). Please refer to the applicable national regulations for the prescribed air replacement rate, max. pressure losses and waste air ducts. The supply and waste air connectors are not prefitted in some models; they are supplied with the cabinet and must be screwed to the bored hole openings in the head part. Please refer to Annex 1 for the position of the supply and waste air connectors.

Models VBF.60.xxx

Bottom cabinets with floor suction (if a base is used) obtain their supply air from beneath the cabinet. If a bottom cabinet without base is used, i.e. is directly positioned on the floor, it will require supply air from the rear wall of the cabinet. In this case, the black rubber stopper must be removed from the supply air inlet.

Models TRG.205.xxx, G90.xxx

In compliance with EN 14470-2 pressurised gas cylinder cabinets **must** be connected to a technical waste air system which leads outside at a safe place.

An air change rate of 10-times is prescribed for combustible and fire-conductive gases and a minimum air change rate of 120-times for toxic gases. The ventilation system must be permanently in operation. The drop in pressure in the cabinet may not exceed 150 Pa. The incoming speed of the air must be at least 20cm/s.

If, however, a cabinet serves the sole purpose of storing inert gases then it needs not be connected to an exhaust air system.

14. Storage volume/combined storage/provision

Please observe the national regulations depending on the type class of the cabinet.

Models TRG.205.xxx, G90.xxx.xxx

The entire volume of the pressurised gas cylinders stored may not exceed 210 l + 10 l for a scavenging gas cylinder.

15. Marking

All models are labelled in compliance with the requirements of EN 14470-1 and EN 14470-2 (in the case of TRG- and G90-models).

16. Opening the cabinet after a fire

Warning: the cabinets may only be opened following a fire after a period of 6 times the duration of the fire. Depending on the duration of the fire, an ignitable vapour-air mixture may have formed. Before opening the cabinets, remove all ignition sources within a radius of at least 10 m around the cabinets. Only open the doors with the utmost care. Only use spark-free tools.

17. Disposal

All models have been constructed so that they may be disposed of in dismantled form, sorted according to materials. The cabinets contain **no** materials which need to be disposed of as special waste.

18. Safety check

Being a safety equipment the cabinets should be checked once a year as minimum for faultless function as per manufacturer's specification.

Please refer to the test plate on the outside of the door for the next inspection date.

This annual inspection can only be performed with the requisite care and to secure your warranty claims in case of fire by an authorised ascos employee (refer also to Splus Agreement in our general catalogue).

We also recommend that you carry out a functional check on a daily and monthly basis.

The daily inspection should also extend to the collecting sumps (any leakage must be mopped up and removed immediately).

The perfect function of the door(s), and the correct fit and state of the fire protection seals should be inspected on a monthly basis. (In case of damage, please contact your specialised dealer to have the cabinet repaired with original parts.)

Many other products for the safe handling of all kinds of hazardous materials can be found in our general catalogue available free of charge by phoning +44 (0) 1952 814 202.

In the case of faults and complaints with regard to our products during and after the warranty period and to request safety inspections please contact your specialised dealer mentioned on the back of your instruction manual.

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